

**HIGH VOLUME SLIDERITE®** 

## **SECTION A**


DETACHMENT


Remove the 4 safety pins by slightly rotating and pulling by the closed end.

Remove the draw latchs by pulling up-out & off of the dewar shell latch catches.

The dewar shell and the Sliderite® pallet can then be detached if required.

Reverse process to reattach.

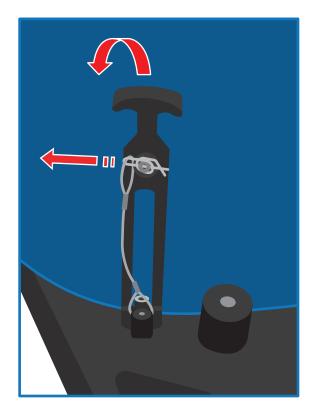

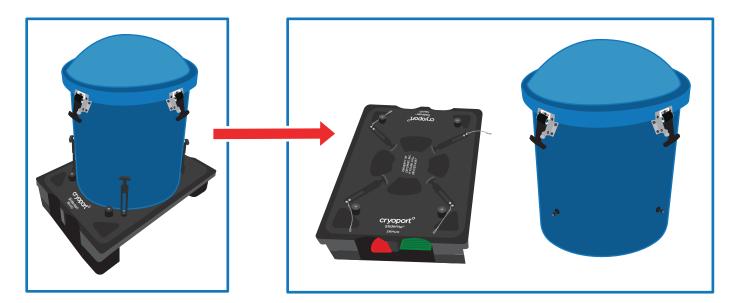

## IF THE DRY SHIPPER DOES NOT EMIT VAPOR WHEN PLUG IS REMOVED PLEASE CONTACT CUSTOMER SERVICE IMMEDIATELY, BY CALLING (949) 470-2305 OR EMAIL CS@CRYOPORT.COM

24-hour Help Line (949) 470-2305 Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614 Phone: 1-949-470-2300 | Fax: 1-949-470-2306 www.cryoport.com • e-mail contact: cs@cryoport.com

ADVANCED THERAPY SHIPPER™