

STEP 1

HOW TO ADD MORE LN2 WITH SAMPLE INSIDE

GENERAL PURPOSE SHIPPER

HIGH VOLUME DRY SHIPPER:


PERSONAL PROTECTIVE EQUIPMENT: CRYOGENIC GLOVES, SAFETY GOGGLES

UNLATCHING INSTRUCTIONS


UNLATCH THE ENCLOSURE PER INSTRUCTION 9356, HOW TO UNLATCH & LATCH HIGH VOLUME ENCLOSURES



Remove the lid and rest it sideways behind the dewar and against the enclosure lid.

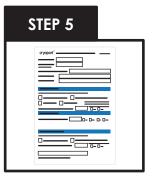

Complete the In Field Recharging form and send it to cs@cryoport.com

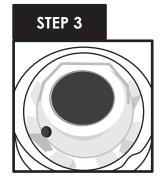

Cryoport recommends the commodity is loaded in a Safepak[®]XL as the bag is capable to withstand LN2. DO NOT REMOVE THE COMMODITY AT ANY TIME.

Slowly pour Liquid Nitrogen into the opening. Cryoport Customer Service will provide information on the amount of LN2 to use for your specific use case.



Replace the lid and let the Dry Shipper stabilize, undisturbed, for at least 30 minutes, Before handing it over to carrier.

IF THE DRY SHIPPER DOES NOT EMIT VAPOR WHEN PLUG IS REMOVED PLEASE CONTACT CUSTOMER SERVICE IMMEDIATELY, BY CALLING (949) 470-2305 OR EMAIL CS@CRYOPORT.COM

24-hour Help Line (949) 470-2305 Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614 Phone: 1-949-470-2300 | Fax: 1-949-470-2306 www.cryoport.com • e-mail contact: cs@cryoport.com