cryoport

SCIENCE, LOGISTICS, CERTAINTY.

Sample Provider

EXP-6 Receiving & Shipping Instructions

As Easy as

Receive Brown Carton



Use caution when using blade to open brown carton. Do not cut through box.





Figure B



Figure C

Remove and discard outer BROWN shipping carton by grasping the top brown carton flap marked with arrows (fig. A) and pulling down along perforations on side of brown carton (fig. B).

Brown carton should be discarded (fig. C). Do not use for any further shipping!

Prepare
White Carton



Flap "A" has paperwork already in place to ship the samples.



LOAD SAMPLES

See "How to Load Samples" instruction sheet for full details.

(Instructions are located on the back side of flap "A")

Ship



When preparing to ship, make sure flap "A" is on TOP of white carton! Be sure you followed "How to Load Samples" instruction sheet.

(Instructions are located on the back side of flap "A")



Seal top of white carton with packaging tape.

Carton is ready for pick-up.

24-HOUR HELP LINE (949) 470-2305 Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614 Phone: 1-949-470-2300 Fax: 1-949-470-2306 www.cryoport.com • E-mail: contact: cs@cryoport.com