

EXP-6 Standard Shipment: Unloading

Step 1



Wear Safety Gloves.
Hold shippo plug in one hand.
Hold blue straps with other hand.
Pull shippo plug upwards and out.

Step 2



Put shippo plug to one side.
Gently pull upwards on blue straps.
Remove sample bag completely
And place on a flat surface.

Step 3



Hold bag with one gloved hand. Use
scissors to cut off end of Tyvek bag.
Remove end of bag completely to re-
veal sample container(s).

Step 4



Remove contents and discard the
Tyvek bag.

Step 5



WARNING! Do not leave any sam-
ples or commodities in the shipper!

Step 6



Replace shippo plug.
Push plug fully into place.

Step 7



If any label was used in this area
for your shipment please place the
enclosed "EMPTY" label over it.

Step 8



Remove Stop Insert from inside shipping envelope on flap "B".
Throw away **ONLY** the Stop Insert.

**Instructions Continued
on Back of Sheet.**

**24-HOUR HELP LINE
(949) 470-2305**

Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614
Phone: 1-949-470-2300 Fax: 1-949-470-2306
www.cryoport.com • E-mail: contact: cs@cryoport.com

Step 9



Close flaps making sure to have flap "B" showing on the outside of the white carton.

FOR INTERNATIONAL SHIPMENTS ONLY:

- REMOVE AND SIGN ALL FOUR (4) COPIES OF COMMERCIAL INVOICES.
- Place all four (4) signed copies of Commercial Invoices back inside shipping envelope on flap "B" behind Air Waybill.

Step 10



Seal top of white carton with packaging tape making sure to have flap "B" showing on the outside of the white carton.

**Please return
empty shipper
to CryoPort
within 24 hours.**

