

GENERAL PURPOSE SHIPPER

EXP6 STANDARD SHIPMENT

PERSONAL PROTECTIVE EQUIPMENT: CRYOGENIC GLOVES, SAFETY GOGGLES



After opening the loaded shipment, ensure the blue Safepak[®] straps stay outside the opening of the dewar at all times.



Wearing safety goggles and cryogenic gloves, pull gently upward to remove the shipper vapor plug.



Cryoport recommends the commodity is loaded in a Safepak[®] as the bag is capable to withstand LN2. **DO NOT REMOVE THE COMMODITY AT ANY TIME.**

STEP 4




- O The amount of LN2 to add to the shipper is provided by Cryoport Customer Service.
- O Slowly pour LN2 into the shipper and avoid spilling on the data logger.
- O The LN2 should absorb into the shipper and excess LN2 should not be visible.

STEP 5





Once the LN2 has been replenished, replace the shipper vapor plug by aligning the slot with the thermocouple wire and gently push the plug into position.

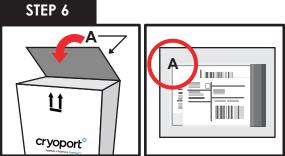

- O Close flaps making sure to have flap "A" showing on the outside of the white carton.
- For International shipments: REMOVE AND O SIGN ALL FOUR (4) COPIES OF COMMERCIAL INVOICES.
- Place all four (4) signed copies of Commercial O Invoices back inside shipping envelope on Flap "A" behind Air Waybill.

STEP 7

Seal top of white carton with packaging tape making sure to have flap "A" showing on the outside of the white carton.



Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614 Phone: 1-949-470-2300 | Fax: 1-949-470-2306 www.cryoport.com • e-mail contact: cs@cryoport.com