

SCIENCE. LOGISTICS. CERTAINTY.

**EXP-6 Shipper** 

# **In-field Recharging Instructions**

# With-out the Commodity Loaded

## PERSONAL PROTECTIVE EQUIPMENT:

- Cryogenic Gloves
  Safety Goggles



## For International or Extended use orders only:

- White sock string could be present, follow Step 2 to discard
- When no string is present, skip Step 2.





Wearing safety goggles and cryogenic gloves, pull directly up o remove shipper plug. Place plug to one side. DO NOT discard plug.

#### Step 2



Pull white string and remove socks.



### **CAUTION!**

There may be residual LN2 in the socks. Hold socks over opening of the Dry Shipper to allow liquid to drip back into the shipper. Use caution when transporting socks to garbage to prevent contact with liquid.

#### Step 3



Slowly pour Liquid Nitrogen into opening. Only pour the amount as instructed by Cryoport Client Care.

#### Step 4

Replace the vapor plug and let the Dry Shipper stabilize, undisturbed, for at least 1 hour.



#### Step 5

Pour off any excess liquid nitrogen by tipping the open end of the dry shipper toward the ground for two seconds.



### Step 6



Replace the shipper vapor plug by aligning the slot with the thermocouple wire and gently push the plug into position.

24-HOUR HELP LINE (949) 470-2305

Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614 Phone: 1-949-470-2300 Fax: 1-949-470-2306 www.cryoport.com • E-mail: contact: cs@cryoport.com

Form: 9200 Ver. 00, Effective Date: 12/02/16