



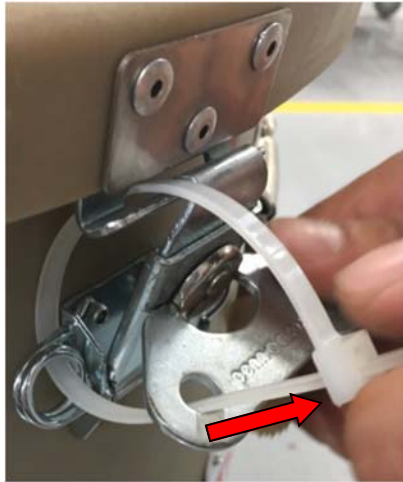
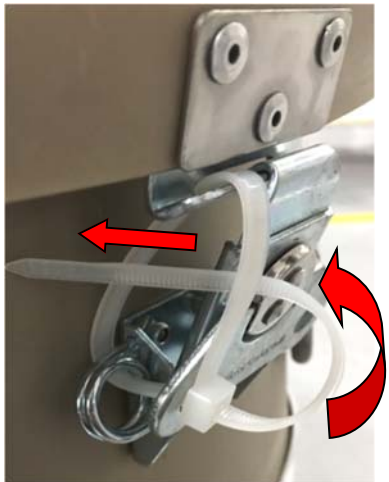


Standard Shipment: CXHVO Latch Zip Tie Procedure

How to Secure CXHVO Latch with Zip Ties

| | | |
|---|---|--|
| <p>Step 1</p> <p>Insert Zip Tie down through Lid Latch Catch</p> |  <p>Assure that it is Through the Latch Catch Slot!</p> | <p>Step 2</p> <p>Thread end of Zip Tie through Body Latch Spring Loop</p>  |
| <p>Step 3</p> <p>Thread Zip Tie Up through Body Latch Turn-buckle hole.</p>  | <p>Step 4</p> <p>Bring the Zip Tie end up to the Buckle</p>  | |
| <p>Step 5</p> <p>Insert Zip Tie through Lid Buckle and tighten</p>  | <p>Step 6</p> <p>Finish by inserting the end of the Zip Tie back through the loop.</p>  | |

24-HOUR HELP LINE
(949) 470-2305

Form: 9210 v1 Effective: 02/20/17

CryoPort, Inc. • 17305 Daimler St. • Irvine, CA 92614
Phone: 1-949-470-2300 Fax: 1-949-470-2306
www.cryoport.com • E-mail: cs@cryoport.com