

SCIENCE. LOGISTICS. CERTAINTY.

Standard Shipment: CXHVO Latch Tamper Proof Tag Procedure for the Shipping Container

Shipping Process


How to Secure CXHVO with Tamper Proof Tag

Step 1

Primary Method:

If utilizing this method install Zip ties to secure Latches Per Instruction 9210




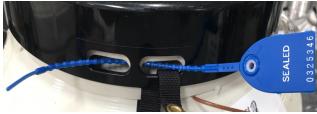



Step 2

Align the slots in the vapor plug lid and the dewar. Close the

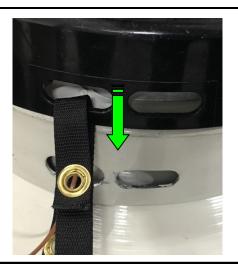

Step 3


Take Tamper resistant Tag ACC-9128 and slightly bend near the tip to aid in inserting the tag though the



Step 4

Insert
Tamper
Tag
through
Vapor
Plug Lid
& Dewar
slots



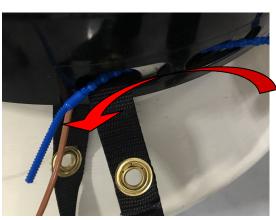

24-HOUR HELP LINE (949) 470-2305 CryoPort, Inc. ● 17305 Daimler St. ● Irvine, CA 92614
Phone: 1-949-470-2300 Fax: 1-949-470-2306
www.cryoport.com ● E-mail: cs@cryoport.com



SCIENCE. LOGISTICS. CERTAINTY.

Standard Shipment: CXHVO Latch Tamper Proof Tag Procedure for the Dewar Body

Shipping Process

How to Secure CXHVO with Tamper Proof Tag

Step 6

Pull the tamper tag though slots until it is centered in the slots.



Step 7


Insert Tamper Tag
Tail
through
Tamper
Tag Flag
retention
hole & Pull
Tight!



Step 1

Alternate
Method: Insert Tamper
Tag down
through Lid
Latch Catch



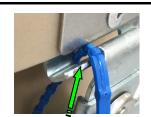

Assure that it is Through the Latch Catch Slot!

Step 2

Thread end of Tamper Tag through Body Latch Spring Loop



Step 3

Thread Tamper Tag Up through Body Latch Turnbuckle hole.



Step 4

Bring the Tamper Tag tip up to the Tamper Tag Flag End



Step 5

Insert Tamper Tag through Tamper Tag Flag retention hole & Pull Tight!



24-HOUR HELP LINE (949) 470-2305 CryoPort, Inc. • 17305 Daimler St. • Irvine, CA 92614
Phone: 1-949-470-2300 Fax: 1-949-470-2306
www.cryoport.com • E-mail: cs@cryoport.com