

**ADVANCED THERAPY SHIPPER™** 

## HIGH VOLUME SHIPMENT: LOADING NON-INFECTIOUS SAMPLES

**SECTION A** 

**RECEIVING SHIPPER** 

#### STEP 1

#### **UNLATCHING INSTRUCTIONS**



Once the shipper has arrived, remove old shipping pouch and zip-ties.



Remove zip-ties from both black handles with scissors.



Unlatch both sides by pulling black handle down & away from the shipper.



Open lid to expose the dewar.

**SECTION B** 

**PACKING THE SPECIMEN** 

IMPORTANT NOTE: Commodity and accessories must be pre-conditioned prior to loading into the sample chamber. A 20% reduction of holding time can occur if the materials are not preconditioned prior to loading into the shipper.

# STEP 1

#### LOADING SAMPLES TO BE SHIPPED INTO EITHER:







**Stainless Steel** Canister 4 canisters can be loaded into the dewar


**Cryovial Box** 

Canes Holds 5 (2ml) vials.



ACC-9139



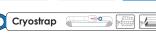

ACC-9141

ACC-9140



Safepak® Holds 15 canes or 2 (5 1/2" x 2" x 2") boxes. 7 bags can be loaded into the dewar.



**Cryostrap®** Use alone or with the Safepak®XL to secure blood cassettes and cryovial boxes





ACC-10502

**Cassette Rack** 

Holds 9 (9.5" x 6 x () 5") cassettes

**Box Rack** 

Holds 5 (5 3/8 x 5 3/8" x 2 1/4") or 10 (5 1/2" x 2" x 2")

ACC-10521

## **SECTION C**

#### LOADING SAMPLES FOR SHIPMENT

# STEP 1



Cut off the zip tie on the hinged cap with scissors.



Open the hinged cap by pulling up on one of the two lift handles on the lid. With gloved hands, pull up on the circular handle in the center of the vapor plug to remove the vapor plug. Set the vapor plug aside.

# STEP 2



Lower the packaged commodity into the dewar with gloved hands.

# STEP 3



If your commodity is loaded in a rack, ensure the handles are laying flat against the top.

#### IF THE DRY SHIPPER DOES NOT EMIT VAPOR WHEN PLUG IS REMOVED PLEASE CONTACT CUSTOMER SERVICE IMMEDIATELY, BY CALLING (949) 470-2305 OR EMAIL CS@CRYOPORT.COM

**ADVANCED THERAPY SHIPPER™** 

#### HIGH VOLUME SHIPMENT: LOADING NON-INFECTIOUS SAMPLES

#### **SECTION D**

#### **PACKAGING THE SHIPMENT**

# STEP 1



Replace the vapor plug.

## STEP 2



Close the hinged cap by pulling down on the lid.

#### STEP 3



Secure the hinged cap with a zip tie.

#### STEP 4



Remove the Leg 2 shipping pouch from the document protector attached to the dewar handle inside of the container.

### STEP 5





Remove any classification labels from the shipping pouch and place on the metal diamond.

#### STEP 6

**Remove the commercial Invoices** located in the shipping pouch.

#### **International Shipments ONLY**



Sign and date.



Return all documents being shipped behind the Air Waybill Including any: Permits, Forms, Licenses, etc.

#### STEP 7

#### Once all documents are loaded:



Close the shipping pouch and remove sticker backing.



Place shipping pouch on the enclosure's metal plate.



Close the enclosure lid.

# STEP 8

#### LATCHING INSTRUCTIONS





Latch handle on both sides by pulling the black handle down & then towards the shipper.



Insert zip-tie through one of the holes on the metal latch hardware.



Thread end of zip-tie through the hole on the other side of the metal hardware.



Insert zip-tie through the lid buckle and tighten.



Zip-tie is now securely around the black handle.

For those shipments containing regulated dangerous goods/hazardous materials, the shipper is responsible for correctly preparing the shipment according to the current International Air Transport Association (IATA) and International Civil Aviation Organization (ICAO) dangerous goods regulations.

This includes correct identification, classification, packaging, dangerous goods markings and labeling as well as completion of all pertinent and required documentation. The shipper is the customer who is preparing the shipment at each leg in the transportation of the dry dewar packaging.

24-hour Help Line (949) 470-2305

Cryoport, Inc. • 17305 Daimler St. • Irvine, CA 92614 Phone: 1-949-470-2300 | Fax: 1-949-470-2306 www.cryoport.com • e-mail contact: cs@cryoport.com