


STANDARD SHIPPER (CXST1)

UNLOADING SAMPLES FROM CRYOPORT EXPRESS® STANDARD SHIPPER (CXST1)

SECTION A

UNLOADING SAMPLES

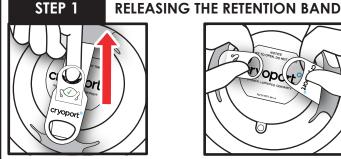

Once the white carton is open, you will need to release the retention band on the shipper plug. Pull the small circle handle up to release the locking tab.



Unlatch and pull apart the straps.



Wear Safety Gloves. Hold the shipper plug in one hand. Hold the blue Safepak[®] straps with other hand. Pull the shipper plug upwards and out.

STEP 3



Put the shipper plug to aside. Gently pull upwards on the blue Safepak® straps. Remove sample bag completely and place on a flat surface.

STEP 4








STEP 5

Remove contents and discard the Safepak[®] bag.



WARNING! Do not leave any samples inside of the shipper. Replace the shipper plug. Push the plug fully into place.

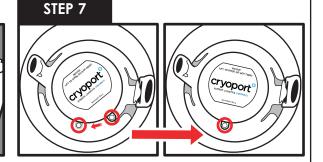

Ensure the hole on the shipper plug and the gray hole on the dewar are aligned. Turn shipper plug to align, if necessary. Also ensure the retention band handles are not overlapping the holes.

STEP 8

container(s).




Use scissors to cut off end of the

completely to reveal sample

Safepak® bag. Remove end of bag

Place the smaller circle handle under the other strap and pull the handle through the slit.

SECURING THE RETENTION BAND

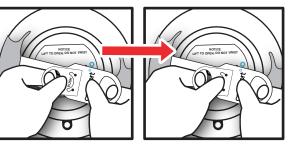

Place your thumb on the word "PUSH" and pull the straps away from each other to place the tab with the word PUSH under the tab with the green check mark.



Ensure the PUSH tab is behind the green check mark tab.



The green check mark will be visible when secured

STEP 9

If you are expecting to receive non-temperature sensitive accessories with your shipment, please see the "How to Include and Remove Accessories in the Standard/Combo Shipper" instructions.

IF THE DRY SHIPPER DOES NOT EMIT VAPOR WHEN PLUG IS REMOVED PLEASE CONTACT CUSTOMER SERVICE IMMEDIATELY, BY CALLING (949) 470-2305 OR EMAIL CS@CRYOPORT.COM

(1/2 fold)




STANDARD SHIPPER (CXST1)

UNLOADING SAMPLES FROM CRYOPORT EXPRESS® STANDARD SHIPPER (CXST1)

SECTION B


PACKAGING THE SHIPMENT

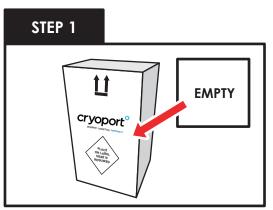

If any label was used in this area for your shipment please place the enclosed "EMPTY" label over it.



Remove Stop Insert from inside shipping envelope on flap "B Throw away ONLY the Stop Insert.

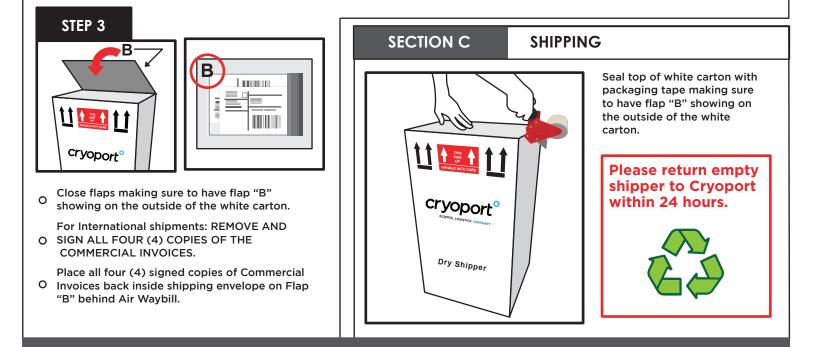

For those shipments containing regulated dangerous goods/hazardous materials, the shipper is responsible for correctly preparing the shipment according to the current International Air Transport Association (IATA) and International Civil Aviation Organization (ICAO) dangerous goods regulations.

This includes correct identification, classification, packaging, dangerous goods markings and labeling as well as completion of all pertinent and required documentation. The shipper is the customer who is preparing the shipment at each leg in the transportation of the dry dewar packaging.

24-hour Help Line (949) 470-2305 Cryoport Systems, LLC • 17305 Daimler St. • Irvine, CA 92614 Phone: 1-949-470-2300 | Fax: 1-949-470-2306 www.cryoport.com • e-mail contact: cs@cryoport.com