


RECEIVING & SHIPPING INSTRUCTIONS

SECTION A



RECEIVE BROWN CARTON



Use caution when using blade to open the brown carton. Do not cut through the box.

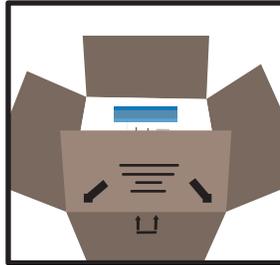

Remove and discard the outer BROWN shipping carton by grasping the top brown carton flap marked with arrows and pulling down along perforations on side of brown carton.

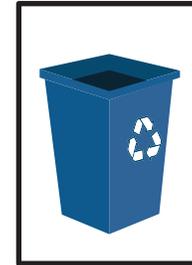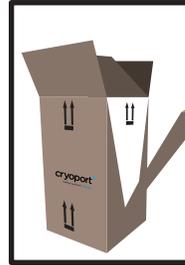

The brown carton should be discarded. Do not use for any further shipping.

SECTION B

PREPARE WHITE CARTON



Flap "A" has paperwork already in place to ship the samples.




See "How to Load Samples" instruction sheet for full details.

(Instructions are located on the back side of flap "A")


SECTION C

SHIP WHITE CARTON

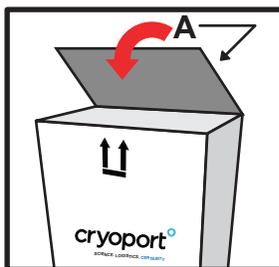

When preparing to ship, make sure flap "A" is on **TOP** of white carton. Be sure you followed "How to Load Samples" instruction sheet.

(Instructions are located on the back side of flap "A")

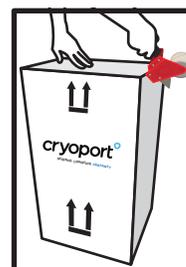

Seal top of the white carton with packaging tape.

Carton is ready for pick-up.

IF THE DRY SHIPPER DOES NOT EMIT VAPOR WHEN PLUG IS REMOVED
PLEASE CONTACT CUSTOMER SERVICE IMMEDIATELY, BY CALLING (949) 470-2305 OR EMAIL CS@CRYOPORT.COM