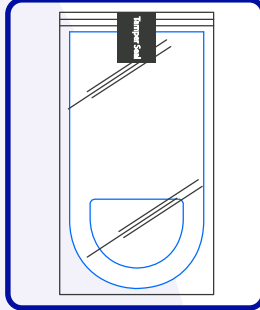


SAFEPAK® XL*Loading and Unloading Instructions*

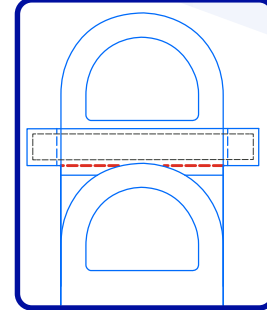
NEED HELP? CONTACT CRYOPORT SYSTEMS' CLIENT CARE AT (949) 470-2305 OR EMAIL CS@CRYOPORT.COM

*Loading Instructions:***Step 1**

Wear the appropriate Personal Protective Equipment (PPE) when handling cryogenic shipments.

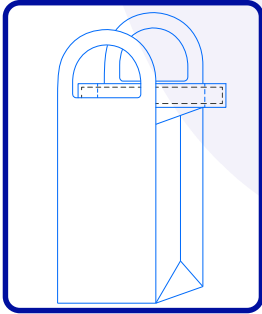
Step 2

Remove the Safepak® XL from the sealed clear bag.

Step 3

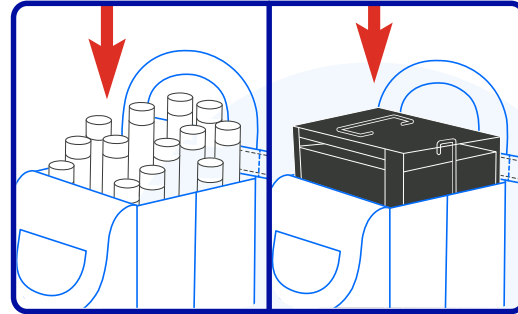
Prefold the top flap along the scored line as highlighted by the **RED** dashed line.

NOTE: The prefold will fold over the adhesive seal strip and behind the handle.

Step 4

To prepare for loading, unfold and open the bag on a flat surface.

NOTE: Fully extend the bag to insert the payload correctly.

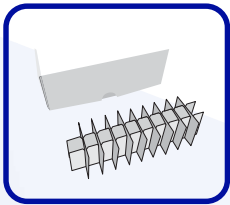
Step 5

Insert cryovial canes, boxes, a loaded metal rack, or other payload into the Safepak® XL bag.

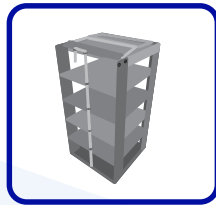
NOTE: Ensure the payload is fully inserted and sitting on the bottom of the Safepak® XL bag to avoid sealing issues.

*Compatible Accessories:***ACC-9043
Cryovial Box**

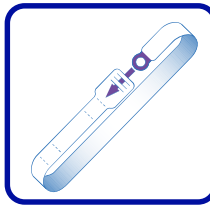
Holds 15 boxes (2" x 5.75" x 2")

**ACC-8905
Box Rack**

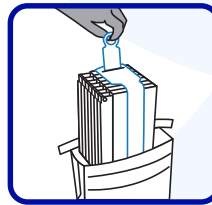
Holds 5 (5.25" x 5.25" x 2") or 10 (5.25" x 2" x 2") boxes

**ACC-9141
Cryostrap®**

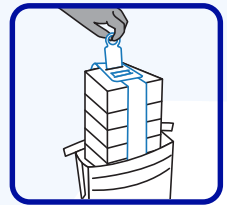
Secures blood cassettes and cryovial boxes

**Cassettes with
Cryostrap®**

Holds 9 blood cassettes (250ml)

**Standard Cryovial
Box with Cryostrap®**

Holds 5 boxes (2" x 5" x 5")



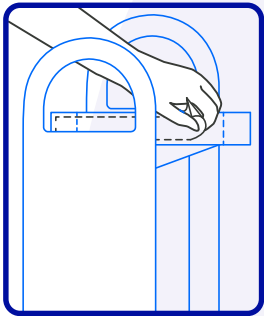
Disclaimer: Safepak® XL can absorb and contain up to 500 ml of liquid.

PROPRIETARY AND CONFIDENTIAL

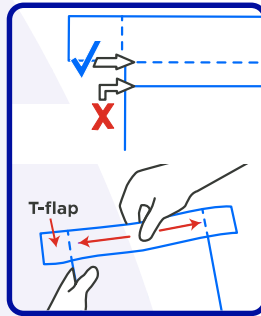
Document retrieved on: 12/26/2024

SAFEPAK® XL*Loading and Unloading Instructions*

NEED HELP? CONTACT CRYOPORT SYSTEMS' CLIENT CARE AT (949) 470-2305 OR EMAIL CS@CRYOPORT.COM

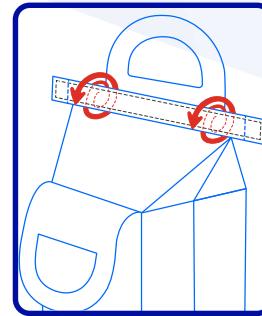
Step 6

Remove the adhesive seal cover from the top flap.

Step 7

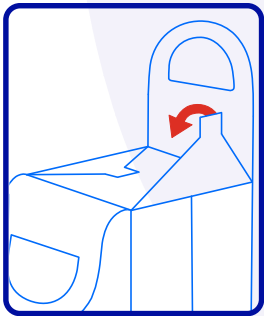
To close the bag, fold down the prefolded score line from Step 3. Then, begin by sealing the **CENTER** of the prefold score line and work your way out to the sides. There should be **NO WRINKLES** or **GAPS** in the seal.

NOTE: Do NOT fold the side T-flaps yet.

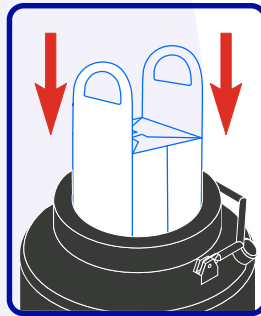
Step 8

Roll the sealed top down the bag until it reaches the top of the payload.

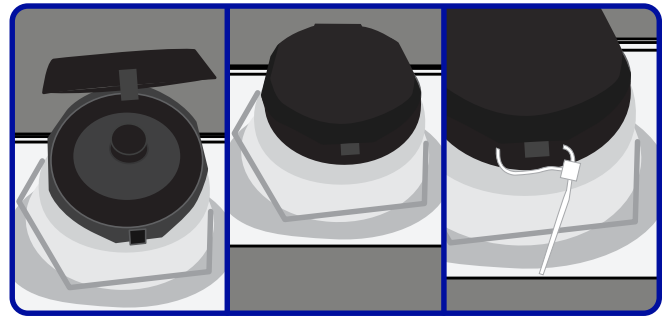
NOTE: Move the handles out of the way when rolling.

Step 9

After rolling the bag down, flatten the sides into a triangle and fold toward the middle. Complete the sealing process by pressing down firmly on the T-flaps, sticking them to the bag.

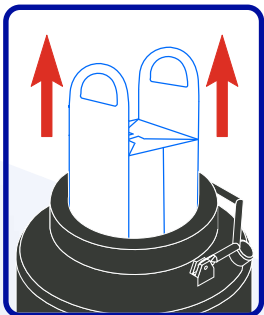
Step 10

Use the handles to lower the packaged Safepak® XL bag into the dewar.

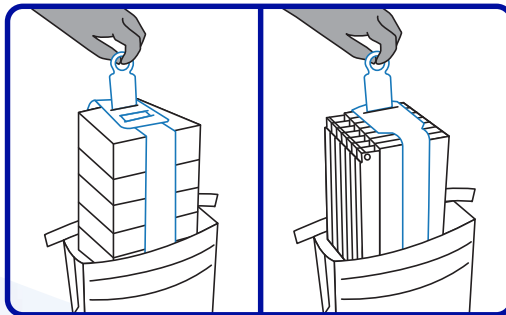
Step 11

Reinsert the vapor plug. Close the hinged cap and secure it with a zip tie.

NOTE: If a serialized zip tie is present, record the serialized number on all necessary documentation.

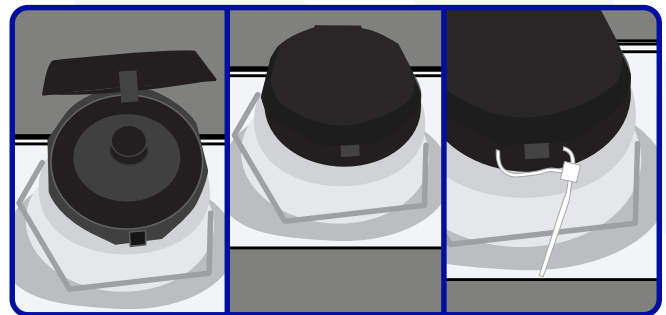
*Unloading Instructions:***Step 1**

Slowly lift the Safepak® XL out of the dewar by pulling upwards.

Step 2

Cut open the top of the Safepak® XL and remove the payload.

NOTE: Please DO NOT return used packaging to Cryoport Systems. Disposal fees will apply accordingly.

Step 3

Reinsert the vapor plug. Close the hinged cap and secure it with a zip tie.

NOTE: If a serialized zip tie is present, record the serialized number on all necessary documentation.

PROPRIETARY AND CONFIDENTIAL